

Data sheet

HigherPurity™ FFPE DNA Isolation Kit

Cat. No: AN0160 (50 reactions)

Description

Formalin fixation and paraffin embedding (FFPE) is a standard method for long-term preservation of most archived pathological specimens. Such samples provide an invaluable resource for subsequent molecular studies of clinical phenotypes, especially genetic studies in which DNA is not available from fresh or frozen tissues.

FFPE tissue is an excellent source of DNA, but its extraction remains a challenge, due to that kind of conservation damage the DNA quality.

Extractions of these types of samples involve two different phases: Deparaffinization and DNA extraction. The procedure omits the use of flammable and malodorous xylene or d-limonene commonly used for deparaffinization, a proprietary buffer formulation DEPARAFFINIZATION SOLUTION is used for the complete dissolution of the wax to release the tissue.

Features

- Silica membrane technology with special MicroSpin columns.
- Very easy paraffin removal.
- Safe method avoids xylene and other toxic.
- Complete removal of contaminants and inhibitors for reliable downstream applications.
- Low elution volume: 15-30 µl.
- The quality of DNA is suitable for the following applications as quantitative PCR or Next generation sequencing (NGS).

Applications

- Rapid isolation of DNA from formalin-fixed, paraffin-embedded samples.
- Isolation of DNA from fresh and archived FFPE samples.
- Isolation of DNA from specimen of object slides.
- Typical downstream application: PCR, qPCR, NGS, STR analysis.

Kit Components

	50 prep
Deparaffinization solution	30 ml
Tissue Lysis Buffer	10 ml
Lysis/ Binding Buffer	15 ml
Proteinase K (*)	2x30 mg
Disinhibition Buffer (**)	18 ml
Wash Buffer (**)	10 ml
Elution Buffer	10 ml
MicroSpin Columns	50 units
Collection tubes	100 units

*Dissolve Proteinase K in nuclease-free water to obtain a 20 mg/mL stock solution(1.5ml/vial). **The Proteinase K solution can be stored for several days at 2–8 °C.** For longer-term storage, the unused portion of the solution may be stored in aliquots at –20 °C until needed.

**Add the volume ethanol (96%-100%) specified [Not included] prior to initial use (see bottle label for volume). After ethanol has been added, mark the bottle to indicate that this step has been completed.

Kit Storage

The kit is shipped at ambient temperature. Upon arrival, store Proteinase K at 4°C, all other kit components can be stored at room temperature. The kit components are stable for 1 year, if stored properly.

Product use limitation

This product is developed, designed and sold exclusively for research purposes. The product was not tested for use in diagnostics or for drug development.

Please refer to www.canvaxbiotech.com for the Material Safety Data Sheet of the product.

(Continued on reverse page)

Distributed by:

Lab Unlimited
CARL STUART GROUP

Tallaght Business Park
Whitestown, Dublin 24,
Ireland
D24 RFK3

Tel: (01) 4523432
Fax: (01) 4523967
E-mail: info@labunlimited.com
Web: www.labunlimited.com

Quatro House, Frimley Road,
Camberley,
United Kingdom
GU16 7ER

Tel: 08452 30 40 30
Fax: 08452 30 50 30
E-mail: info@labunlimited.co.uk
Web: www.labunlimited.co.uk



DETAILED PROTOCOL



A) DEPARAFFINIZATION SAMPLE

1. Add **400 µl of Deparaffinization solution** to one FFPE tissue slice (10µm). **!! Make sure the sample does not comprise more than 15 mg paraffin. Vortex for 10 seconds.**
2. Incubate **3 minutes at 60°C** to promote the melting of the paraffin.
3. Vortex the sample immediately (at 60°C) at a vigorous speed to dissolve the paraffin. After 3 minutes you will see the tissue floating in the wax remover.
4. **Centrifuge full speed for 3 minutes** to pellet the tissue.
5. **Remove the supernatant** with pipetting, avoiding the tissue.
6. **Add 1 ml ethanol 100%**. Vortex for 20 seconds.
7. **Centrifuge full speed for 3 minutes** to pellet the tissue. Remove the ethanol with pipetting, avoiding the tissue.
8. Place the tubes **at 55°C for 10 minutes** with caps open to evaporate the ethanol.

B) DNA ISOLATION

9. Add **200 µl of the Tissue Lysis Buffer + 40µl Proteinase K**. Mix by vortexing 2-5 seconds. **Incubate at 55°C** with agitation (600 rpm) overnight or until the tissue is completely dissolved.
10. Add **20µl Proteinase K** and incubate for 1–2 h at 55°C. Centrifuge at 8000 rpm for 60 seconds.
11. **Transfer the supernatant** into new 1.5-mL microcentrifuge tube.
12. **Incubate at 90°C for 1 hour.** *This incubation partially reverses formaldehyde modification of nucleic acids. If using only one heating block, leave the sample at room temperature after the 55°C incubation until the heating block has reached 90°C.*
13. **Briefly centrifuge** to remove drops from the inside of the lid.
14. Add **300µl of Lysis/Binding Buffer**. Mix by vortexing. **Incubate at Room temperature for 5-10 minutes.**
15. Add **100µl of Isopropanol to the lysate**. Mix by vortexing. Briefly centrifuge the 1.5 ml tube to remove drops from inside the lid.
16. **Transfer the lysate** into reservoir of a combined MicroSpin Column –collection tube.
17. **Centrifuge at 8000 rpm for 60 seconds.** Remove the collection tube. If the sample is not drawn completely through the matrix, repeat the centrifugation step.
18. Place the MicroSpin column in a new collection tube and add **500 µl of Disinhibition Buffer.**
19. **Centrifuge at 12000-14000 rpm for 60 seconds.** Remove the liquid.
20. **Add 500 µl of Wash Buffer** into reservoir of MicroSpin column.
21. **Centrifuge at 12000-14000 rpm for 60 seconds.** Remove the liquid.
22. **Centrifuge at maximum speed for 3 minutes to remove the residual ethanol.**
23. Remove the collection tube and insert the MicroSpin column in a 1.5 ml microtube. **Add 15- 30µl of Elution Buffer** (preheated at 70°C) directly onto the centre of the silica membrane.
24. **Incubate 1 minute. Centrifuge at maximum speed for 60 seconds.** The microtube contains now genomic DNA.

This product is developed, designed and sold exclusively for research purposes.

Distributed by:

Lab Unlimited
CARL STUART GROUP

Tallaght Business Park
Whitestown, Dublin 24,
Ireland
D24 RFK3

Tel: (01) 4523432
Fax: (01) 4523967
E-mail: info@labunlimited.com
Web: www.labunlimited.com

Quatro House, Frimley Road,
Camberley,
United Kingdom
GU16 7ER

Tel: 08452 30 40 30
Fax: 08452 30 50 30
E-mail: info@labunlimited.co.uk
Web: www.labunlimited.co.uk