

Retsch

Solutions in Milling & Sieving

# We assist you in your daily work

The "Assistants" form the perfect complement to the range of mills and sieve shakers from RETSCH. They optimize your results and increase your working efficiency – benefits that you should take advantage of.

## **Dividing** – as important as the analysis itself!

Errors in sample preparation, above all in dividing the laboratory sample, accumulate in the course of the process and cannot be compensated, even by the most upto-date analytical instruments.

# Only the exact division of the entire sample into representative fractions will guarantee exact and unbiased analytical results!

For quality-conscious manufacturers, processors and supervisors, this means that an exact sample division must be as important as an error-free analysis. Just how large the error can be with different sampling and sample division methods is shown in the diagram below.

RETSCH offers sample dividers which operate with the most exact division methods:

■ Sample Divider PT 100 for dividing the whole sample into representative individual fractions

More on page 4

Rotary Tube Sample Divider PK 1000 for extracting representative individual samples from a material flow

More on page 6

 Sample Splitter RT for economically priced manual on-site sample dividing More on page 7

## Important rules for sample division:

#### 1. Significance of the minimum divided quantity

For the sub-sample to be representative of the entire bulk, its qualitative properties must correspond to that of the original quantity. In order to comply with this requirement, some basic guidelines should be respected. In particular, this concerns the minimum partial quantity as well as the maximum permissable deviation. There are several application-related guidelines and specifications as to the calculation of the minimum quantity and sampling accuracy. Please refer to DIN 51 701, part 2, or the "Guidelines of Preparing Laboratory Samples" from AAFCO (Association of American Feed Control Officials Inc.).

#### 2. Observance of the sample size opening

With all sample dividers used, the sample size opening should be approx. 2.5 – 3 times larger than the maximum particle size. This ensures that even the larger particles can pass through the sample openings uniformly.


#### 3. Exclusion of manual influences

Care must be taken that neither the division process nor the division results are altered by manual influences. This means that the entire fraction obtained must be used for further processing or subsequent analyses!

# Comparison of different sampling and sample division methods

The diagram shows how large the error can be for different sampling and sample division methods. It can clearly be seen that rotary tube sample dividers produce the smallest qualitative variation (A). They achieve the highest degree of fractionating accuracy and are therefore clearly superior to all other methods. Sample splitters provide the best results of all the manual fractionating methods.

- A: Rotary tube sample dividers
- B: Sample splitters
- C: Disc dividers
- D: Heaping and dividing
- E: Random sampling (e.g. with a scoop)

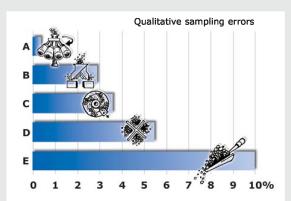

Example: Bulk material, feed size < 5 mm

## Feeding - reproducible and efficient

The controlled continuous material feed with the Vibratory Feeder DR 100 offers many advantages. Errors resulting from discontinuous manual feeding are avoided and the results are always reproducible. The DR 100 can be used for material feeding in combination with sample dividers, mills and many other instruments.



The benefits of automatic feeding:

- Uniform feed of larger amounts
- No temporary overloads
- Less time- and labor-intensive
- Reproducible results

The Vibratory Feeder DR 100 can be used in many applications. The wide range of accessories makes the instrument adaptable to your own particular requirements.

More on page 9

**Drying** – rapid and gentle

The drying of materials in the RETSCH Fluid Bed Dryer TG 200 takes place in a constant flow of warm air. The result is a rapid, gentle and uniform drying process. The savings in time compared with drying in a drying oven are considerable and may be up to several hours. After drying in the fluidized bed the bulk goods are both pourable and thoroughly mixed.



Many substances have already been successfully dried in the TG 200, e.q.:

- Chalk, coal, coke, compost, fertilizers, grain, hops, ion exchangers, metal powders, peat, plant parts, pharmaceutical products, plastics, polymers, rubber, salts, sand, sawdust, soil, tea, tobacco, washing powder, waste.
- Apart from bulk goods it is also possible to rapidly dry test sieves which can be installed directly on the instrument.
   More on page 10

## Cleaning - care is required



RETSCH's product range includes three sizes of ultrasonic baths for cleaning test sieves and grinding tools quickly and easily. UR 1 is for test sieves up to 203 mm dia., UR 2 for test sieves up to 400 mm dia., and the UR 3 for the simultaneous cleaning of up to 5 test sieves 200/203 mm dia. The gentle yet thorough cleaning of test sieves in an ultrasonic bath

increases their working lives as damage which could occur during manual cleaning is avoided.

More on page 12

Do you need any further "Assistants"?

Please let us have your proposals about how we can make working with RETSCH mills and sieve shakers even more convenient for you. We will be grateful for your suggestions!

# Pelletizing - Sample preparation for XRF

For the preparation of solid samples for XRF analysis RETSCH offers pellet presses in 2 sizes. The automatic floor model PP 40 features individual pressure force regulation up to 40 t. The pellets are pressed into steel rings with external diameters of 40 or 51.5 mm. A version with aluminum cups is also available. The manual hydraulic Pellet Press PP 25 is a compact benchtop unit with pressing tools for 32 mm and 40 mm pellets.

More on page 14



# Sample Divider PT 100



# Simply representative division results

RETSCH sample dividers are rotating dividers. They divide all pourable solids up to 10 mm so accurately that the characteristic composition of each fraction of the sample corresponds exactly to that of the original bulk sample. This very high degree of dividing accuracy and reproducibility is achieved with both fine and coarse materials. The material feed and dividing processes take place automatically, without interruption and without loss of material. The feed amount can range from a few grams up to 5000 ml depending on the sample vessels used. It is possible to produce an individual number of identical fractions for various applications by the repeated division or combination of fractions.

## Benefits at a glance

- Extremely high division accuracy
- Representative and reproducible division for accurate analysis results
- Modular design
- Automatic material feed via synchronized feeder
- Simple and rapid handling due to a convenient quick-release clamping system for sample vessels
- Speed is monitored and kept constant
- Timer
- Compact, maintenance-free and easy to clean
- Low-noise drive
- CE-conforming
- 2-year warranty

# The easy way of sample division

Working with the RETSCH Sample Divider PT 100 is easy and uncomplicated. For example, material feed with the Feeder DR 100 is automatic and synchronized: this means representative sample division right from the start, as the DR 100 only starts feeding the material when the sample divider has reached its proper running speed.

For cleaning and exchanging, the dividing head, feed chute and hopper can be easily removed without any tools. The sample vessels are also extremely easy to attach and release with the quick-release clamps. This new type of clamping system no longer involves tiresome mounting of the vessels.

10013.	
Performance data	PT 100
Applications	sample division, sample reduction
Feed material	bulk materials
Number of divisions	6, 8 or 10
Timer	1 - 60 min or continuous
Technical data	
Feed size	≤10 mm
Feed canacity	may 5000 ml

 Feed size
 ≤10 mm

 Feed capacity
 max. 5000 ml

 Vessel volume
 30, 100, 250 or 500 ml

 W x H x D
 580 x 910 x 420 mm (incl. DR 100)

 Net weight
 approx. 33.5 kg (incl. DR 100)

Noise values (Noise measurement according to DIN 45635-31-01-KL3)

Measuring conditions: Dividing material: silica sand; particle size <3 mm Emission value with regard to workplace  $L_{pAeq} \ 41 \ dB(A)$ 

### Versatile, variable, individual – sample dividing with the RETSCH PT 100

The Sample Divider PT 100 has a modular design and can be put together to suit individual requirements. It offers an extremely flexible range of possible applications. A feeder, various dividing heads, sample receptacle vessels and further useful accessories are available in addition to the drive unit.

The number of part samples is deter-

mined by the choice of the dividing

Sample vessels are available in different sizes for various applications. Wide-mouth glass bottles fit the dividing heads as standard. Special dividing heads are available for use with Duran laboratory bottles (100, 250 and 500 ml). These dividing heads can also be equipped with inserts for 30 ml plastic bottles.

the use of a protective cap for the dividing head hopper. The dust cap minimizes both material losses and dust formation.

For fractions with a low density or with a high fineness we recommend

For uniform material feed the Vibratory Feeder DR 100 should be used. PT 100 and DR 100 are connected via an interface and therefore perfectly matched. In addition, the special swivel back stand assures a fixed position of the feeder over the center of the dividing head inlet thus ensuring increased accuracy. Further information about the Feeder DR 100 can be found in this brochure on page 9.

RETSCH offers a complete unit which includes an 8-outlet aluminum dividing head with the practical quick-release system. The set is supplied with 10 wide-mouth 250 ml sample bottles.

Order data on page 8



- Dividing head with quick-release system for wide-mouth bottles
- Dividing head with quick-release system for particle sizes <5 mm, for use with Duran bottles and for
- Insert for 30 ml plastic beakers

# **Technology of PT 100**

The material to be divided first flows through a decentrally located feed hopper directly into the openings in the dividing head. Even with coarse material, this achieves a very low level of deviations between the materials in the sample vessels. The dividing process itself runs automatically and without manipulation. The dividing head rotates – with speed monitoring – at a constant 110 revolutions per minute, independently of the load and the

mains frequency. That means that with a dividing head with ten outlets, the feed flow is divided into 1100 individual samples each minute. The highest degree of dividing accuracy is thus guaranteed. The dividing heads divide the material evenly among the sample vessels. Depending on the quantity and further application, amongst others wide-mouth bottles and Duran laboratory bottles can be used.



# **Rotating Tube Divider PK 1000**



# Benefits at a glance

- Exact dividing, also of larger quantities
- Representative and reproducible division for accurate analysis results
- Modular design
- Adjustable dividing ratio
- Extraction of 1-3 samples
- Dividing process according to DIN 51 701/Pt 4
- Batch and continuous operation possible
- Easy cleaning
- 2-year warranty

# Easily divides large quantities

The RETSCH rotating tube divider is the prerequisite for representative dust-free division and volume reduction of larger bulk samples. It is suitable for powdered or granular bulk materials with particle sizes up to 10 mm. The rotating tube divider can be provided with bottom cones for 1, 2 or 3 samples. The slot width adjusts the ratio of the fractions and therefore the amount of sample. The minimum feed amount should not be less than 100 ml in order to ensure maximum accuracy.

The sample fractions can be collected in laboratory bottles with a capacity of up to 0.5 liters. The reject collector has a capacity of 30 liters.

All parts coming into contact with the sample material are made from stainless steel or glass. In order to ensure uniform material feed and therefore an increased sample division accuracy, we recommend that the rotating tube divider be used in conjunction with a DR 100 feeder.

The dividers are also suitable for inclusion in continuously working laboratory and pilot-plant scale units.

The PK 1000 is available as a complete unit including 10 x 500 ml sample bottles, a 30 l reject collector, bottom cone and vibratory feeder. It is also possible to select the components individually according to your particular requirements.

Performance data		PK 1000	
Applications	samp	ling and sample di	viding
Feed material		bulk materials	
Number of divided samples		1 - 3	
Timer	digital,	1 - 99 min or con	tinuous
Technical data			
Available bottom cones	with 1 sample outlet	with 2 sample outlets	with 3 sample outlets
Slot width, continuously adjustable	0 - 159 mm	0 - 110 mm	0 - 53 mm
Max. dividing ratio	1 x 1:5	2 x 1:7.2	3 x 1:15
Min. dividing ratio*	1 x 1:26	2 x 1:26	3 x 1:26
Feed size	≤10 mm	≤10 mm	≤10 mm
Volume of reject collector		30 liters	
WxHxD	56	60 x 1150 x 700 m	ım
Net weight	approx	. 37 kg (without D	R 100)

\* for a maximum particle size of 10 mm. for smaller maximum particle sizes the division ratio increases accordingly

Noise values (Noise measurement according to DIN 45635-31-01-KL3)

Measuring conditions: Dividing material: silica sand; particle size <3 mm Emission value with regard to workplace  $L_{\text{DAe}\alpha} \ 63 \ dB(A)$ 

#### **Technology of PK 1000**

The material to be divided passes through the feed hopper into the rotating tube divider. The total material stream is distributed evenly at a constant speed (50 min<sup>-1</sup>) over the pitch circumference of the lower cone by the tube rotating within the upper cone. The interchangeable lower cones have one, two or three continuously adjustable sample

slots. In the course of each rotation a separated quantity corresponding to the width of the slot is deposited in the sample bottle. The rest passes into the reject collector.



### Calculating the slot width for the PK 1000

The maximum dividing ratio depends on the maximum slot width which can be set on the bottom cone. This differs between the three bottom cones which are available (see table). The smallest dividing ratio depends upon the maximum particle size of the sample since the slot width should be at least 3 times wider than the maximum particle size.

This means that smaller fractions can be taken from smaller particle sizes. The slot width "X" to be set can be calculated from the ratio of the required fractional amount "OT" to the initial sample amount "QA" multiplied by the fixed pitch circumference "U" of the bottom cone (U = 795 mm for all bottom cones).

$$X = \frac{QT}{QA} * U$$

Example: A representative sample of 250 ml is to be taken from an initial sample amount of 5000 ml. This means that the slot width must be set to 40 mm.

Order data on page 8



# Sample Splitters RT 6.5 - RT 75

# **Accurate manual dividing**

RETSCH sample splitters are used for the simple dividing and reduction of bulk materials of all kinds. Sample



splitters are ideal for the on-site reduction of sample material. They are easy to use, easy to clean and do not need an electrical power supply. Depending on the particle size, material and particle size distribution, the opening width of the passage should be 2.5 - 3 times greater than the diameter of the largest particle (particle size factor). Each sample splitter consists of one dividing head, one stand and three receptacles.

#### Order data on page 8

# Benefits at a glance

- For use in the laboratory and on-site
- High-precision manual dividing process
- Easy and quick to clean
- Dividing process according to DIN 51701, Part 4
- Inexpensive
- Available in 6 sizes

Technical data	RT 6.5	RT 12.5	RT 25	RT 37.5	RT 50	RT 75	
Slot size	6.3 mm	12.5 mm	25.0 mm	37.5 mm	50 mm	75 mm	
Number of slots	12	18	16	12	8	6	
Max. feed size*	approx. 4 mm	approx. 8 mm	approx. 16 mm	approx. 25 mm	approx. 33 mm	approx. 50 mm	
Max. feed charge	5 lit	ters		16 liters			
Material of dividing head	stainles	ss steel	sheet steel, hot-dip galvanized				
Material of stand	sheet stee	el, painted	sheet steel, hot-dip galvanized				
Material of receptacles	tin p	tin plate		sheet steel, hot-dip galvanized			
WxHxD	300 x 270	x 250 mm	620 x 420 x 260 mm				
Net weight	approx	. 3.5 kg		approx.	21.5 kg		

st with a 5-10% fraction of the maximum particle size

#### **Technology of sample splitters**

With sample splitters, the sample material is evenly distributed in one of the receptacles and then emptied over the dividing head. The material runs through the alternately arranged passages in opposite directions into the two collecting receptacles under the dividing head outlets. With every

operation the feed sample is split in halves. This can be repeated as many times as necessary until the required dividing quantity has been obtained.



# Order data sample dividers

Sample divider PT 100						Item No.
PT 100 complete unit, incl. dividing head with 8 quick-release outlets, DR 100/40* and 10 wide-mouth bottles 250 ml						
PT 100 complete unit for 220-240 V, 50 Hz						40.534.0005
PT 100 complete unit for 110-120 V, 6	50 Hz					40.534.0006
PT 100 drive unit (without dividing he	ad, vibratory feede	r/stand, sample recep	otacles)			
PT 100 drive unit for 100-120/200-24	0 V, 50-60 Hz					40.534.0002
Dividing heads for PT 100						
material	sample outlets	feed size	sample bottle		bottle fixation	
aluminum hard-anodized	6	up to 10 mm	wide-mouth bott	les	quick-release	42.793.0003
aluminum hard-anodized	8	up to 10 mm	wide-mouth bott	les	quick-release	42.793.0001
aluminum hard-anodized	10	up to 10 mm	wide-mouth bott	les	quick-release	42.793.0002
POM	8	up to 10 mm	wide-mouth bott	les	quick-release	42.793.0007
aluminum hard-anodized	8	up to 5 mm	Duran laboratory	bottles	quick-release	42.793.0009
Support with 30 ml plasti	c beakers and lids,	8 pieces (for dividing	head 42.793.000	9)		42.018.0001
Sample vessels for PT 100			30 ml	100 ml	250 ml	500 ml
Wide-mouth bottles, 10 pieces			-	-	22.523.000	22.523.0002
Duran laboratory bottles, 10 pieces			-	22.523.000	3 22.523.000	22.523.0005
Plastic beaker with cover 30 ml, 10 pi	eces (for support 4	2.018.0001)	42.156.0001	-	-	-
Accessories for PT 100						
Hopper, stainless steel, volume 2.8 l (	for PT 100 used wi	thout DR 100)				03.785.0146
Protective dust cap for hopper, of PON	1					03.742.0013
Stand for Vibratory feeder DR 100 on	PT 100					42.742.0011

Rotating Tube Divider PK 1	Item No.			
PK 1000 complete unit, incl. bottom c				
10 wide-mout	h bottles 500 ml and reject col	lector 30 liters		
PK 1000 complete unit for 220-240 V,	50 Hz			40.411.0010
PK 1000 complete unit for 100-120 V,	60 Hz			40.411.0011
PK 1000 drive unit (without bottom co	one and vibratory feeder, with 1	0 wide-mouth bottles 500	ml and with reject collector	r 30 liters)
PK 1000 drive unit for 220-240 V, 50/	60 Hz			40.411.0001
PK 1000 drive unit for 100-120 V, 50/	60 Hz			40.411.0006
Bottom cone for PK 1000	sample outlets	slot width	max. dividing ratio	
	1, adjustable	159 mm	1: 5	42.787.0001
	2, adjustable	110 mm	2 x 1: 7.2	42.787.0003
	3, adjustable	53 mm	3 x 1:15	42.787.0004
Sample vessels for PK 1000				
Wide-mouth bottles 250 ml, 10 pieces	3			22.523.0001
Wide-mouth bottles 500 ml, 10 pieces	\$			22.523.0002
Accessories for PK 1000				
Spare reject collector, 30 l				05.010.0004

Sample splitters RT 6.5 - RT 75	Item No.
Sample splitters RT 6.5 and RT 12.5 (incl. 3 receptacles 2.5 liters each, stand and dividing head)	
Sample splitter RT 6.5 with 12 slots 6.3 mm	40.610.0001
Sample splitter RT 12.5 with 18 slots 12.5 mm	40.610.0002
Sample splitters RT 25, RT 37.5, RT 50 and RT 75 (incl. 3 receptacles 8 liters each, stand and dividing head)	
Sample splitter RT 25 with 16 slots 25.0 mm	40.610.0003
Sample splitter RT 37.5 with 12 slots 37.5 mm	40.610.0004
Sample splitter RT 50 with 8 slots 50.0 mm	40.610.0005
Sample splitter RT 75 with 6 slots 75.0 mm	40.610.0006
Spare parts for sample splitters	
Spare receptacle 2.5 liters (for RT 6.5 and RT 12.5)	42.147.0001
Spare receptacle 8.0 liters (for RT 25, RT 37.5, RT 50 and RT 75)	42.147.0002

<sup>\*</sup>Accessories for DR 100 see on page 13

# Vibratory Feeder DR 100

#### Uniform, continuous feeding

The RETSCH vibratory feeder is used for the uniform, continuous feeding and conveyance of pourable bulk materials and fine powders. The DR 100 feeds RETSCH mills and sample dividers, as well as balances and particle measuring devices, and it is also suitable for filling and dosing. Their performance, adaptability and compact design make these devices suitable for a great variety of applications. The DR 100 can also be driven and controlled externally via the built-in interface.

RETSCH vibratory feeders guarantee reproducibly exact results and the economic use of downstream laboratory and testing devices. The DR 100 is easy to set up and operate.

The RETSCH vibratory feeders are available in various designs. For the feed of pourable powders and fine grained bulk materials, we recommend feeding kits with 15 mm chute width, and for granulates and coarser materials feeding kits with 40 mm or 75 mm chute width. The conveying rate is continuously adjustable. Depending on the feed material, device and setting, it can amount to approx. 5 l/min.

For special applications coated chutes, a vibrating tube and a V-shaped chute are available. Aluminum chutes are particularly suitable for samples containing fat or oil.

Order data on page 13

Examples of use
1. DR 100 with Sample Divider PT 100
2. DR 100 with Sample Divider PK 100
3. DR 100 with Ultra Centrifugal Mill ZM 200
4. DR 100 with CAMSIZER from RETSCH Technology









# Benefits at a glance

- Uniform material feed for reproducibly exact results
- Infinitely adjustable volume flow
- Material bed level can be variably adjusted
- Optional control via external interface
- Digital time setting
- Compact control and feed unit
- Easy cleaning of push-fit feed chute
- Maintenance-free design to CE standard
- 2-year warranty

	- T W				
Performance data			DR 100		
Applications		feeding, conveying			
Feed material		pou	rable bulk mate	rials	
Time display		digital, 1	l - 99 min or co	ntinuous	
Volume flow		0 - 5 l/r	min, infinitely ad	justable	
Technical data					
Available chute widths		15 mm	40 mm	75 mm	
Chute length		210 mm	210 mm	210 mm	
Hopper volume		2.8 liters	2.8 liters	3.5 liters	
Feed size		<2 mm	<6 mm	<12 mm	
WxHxD		26	0 x 420 x 280 n	nm	
Net weight			approx. 10 kg		
Noise values (Noise	: measurement a	ccording to DIN	I 45635-31-01-	KL3)	
Measuring conditions: Co	nveying material:	silica sand; part	icle size <1 mm	1	

#### **Technology of DR 100**

The feed material passes through the hopper onto the vibrating chute. This is made to vibrate at 50 (or 60) Hz by two electromagnets. The volume flow can be continuously adjusted. Via the hopper adjustment, the layer height can be set as required. The DR 100 can be driven externally via an interface for example, when it is used in

combination with the Ultra Centrifugal Mill ZM 200. Then the volume flow is automatically adapted to the grinding capacity of the mill. Due to their compact, maintenancefree design the vibratory feeders can be integrated simply into many devices or laboratory installations.

Emission value with regard to workplace



LpAeq 36 dB(A)

# Fluid Bed Dryer TG 200



#### Benefits at a glance

- Gentle drying, dispersing and mixing also of temperaturesensitive materials
- Very short drying times
- Digital parameter setting
- 9 parameter combinations can be stored
- Interval operation
- Versatile with a choice of drying containers and exhaust air filters
- Easy handling with clamping device "comfort"
- Motor without brushes allows for long service life
- Conforming to CE requirements
- 2-year warranty

## Rapid and gentle fluidized bed drying in the laboratory

The dryer TG 200 is used in quality control, sample preparation and R&D departments. It permits the gentle drying of organic, inorganic, chemical or pharmaceutical bulk materials **without localized overheating**. Suitable materials can be coarse, fine, crystalline, fibrous or leafy. The powerful fan ensures optimal air throughput so that the products to be dried are loosened up and thoroughly mixed. With the interval operation the fluidized bed is mixed even better. Temperature, drying time and air volume can be set digitally and adjusted continuously.

In comparison to conventional drying ovens or microwaves, the fluidized bed drying of the TG 200 results in a considerably better performance. The fan produces an air volume of 185 m³/h in idle speed. **The average drying time lies between 5 and 20 minutes**, depending on the type, amount and moisture content of the material. This represents a substantial saving in time and is also favorable for the product which is exposed to less thermal stress.

# The TG 200 is suitable for the following applications:

- Drying of sample materials such as coal, fertilizer, plant parts, plastics, recycling wood, sawdust, secondary fuels, soils and waste. As the motor is located outside the filtered air flow, the TG 200 can also be used for drying more sensitive materials, like e.g. pharmaceutical products, without the risk of sample contamination
- Drying of test sieves

The delivery scope of the fluid bed dryer includes a clamping device "comfort" with filter bag. It is used for attaching the 6 l drying container. Test sieves with 200 mm diameter are mounted directly on the TG 200 without using the drying container. (Adapter for 8" /203 mm sieves available on request).

Drying test sieves with the TG 200

Performance data	TG 200			
Applications	drying			
Feed material	bulk materials and solids, >63 μm			
Temperature control	continuously adj., 40 - 150 °C (dependent on air throughput rate)			
Time setting	continuously adjustable, 0 - 99 min, continuous operation			
Drying time	5 - 20 min, depending on product, quantity, moisture content			
Container volume	1 x 6 liters or 3 x 0.3 liters			
Technical data	TG 200			
WxHxD	400 x up to 1000 x 480 mm			
Net weight	approx. 21 kg			
Noise values (Noise measurement according to DIN 45635-31-01-KL3)				
Measuring conditions: Dried	l product: clay; max. heating power; max. air volume			
Emission value with regard	to workplace L <sub>DAeq</sub> 75 dB(A)			

#### Accessories for the TG 200

6 I drying container, glass or stainless steel, with base made from Conidur stainless steel perforated plate with 63 μm holes. The glass container makes the degree of dispersion of the material during the drying process visible. Thus, the operator can directly adjust the air flow if necessary.

# Clamping device "comfort" with replaceable filter fleece insert

The filter fleece is mainly used for samples with a particle size below  $100~\mu m$ . It allows for sample recovery with minimal loss. The filter can be quickly and easily replaced after each application.



Clamping cover with replaceable filter fleece insert



#### Attachment with 3 removable glass containers (each 0.3 I)

It permits the simultaneous drying of three samples, also of different materials, under the same conditions. This helps to avoid cross contamination. The glass containers can be easily locked and released with a single turn. The perforated plate is made from stainless steel. Container lids with filter fleece inserts are available as optional accessories.

Order data on page 13

## **Technology of TG 200**

Drying in the fluid bed dryer makes use of the fluidized bed process, a technique similar to the one used in large industrial dryers. Ambient air is drawn in through a filter. A blower moves the air across the heating elements, and ultimately forces it through the perforated plate and into the detachable drying container. The solid particles are blown upward and agitated and thus kept separate one from another. This helps to avoid a caking and sticking of the particles as it often occurs when other drying methods are used.

The air stream extracts moisture from the particles and then exits through the filter bag in the cover. Using the quick-clamp cover with the filter fleece insert is advisable when dealing with products finer than 100 µm in diameter.

The 1000 watt blower provides an air volume of 185 m³/h at idle speed; heater output is 2000 watts. The air volume, heating power and temperature are infinitely adjustable. Temperature control is effected using the display gauge.



# Ultrasonic Baths UR 1 / UR 2 / UR 3



### Cleaning

UR 1

The RETSCH ultrasonic baths clean test sieves, microprecision sieves, glass and metal components as well as metallurgical and geological samples, spectacles, jewellery or coins gently and intensively. In addition to cleaning, the ultrasonic baths can also be used for other working processes.

# Benefits at a glance

- Fast, gentle, and highly efficient cleaning
- Universal and compact
- Easy to use, saves time and cost
- Intensive dispersion and degassing
- Low-maintenance, long life and environmentally sound
- 2-year warranty

### **Dispersion**

The RETSCH ultrasonic baths are used to prepare suspension samples for wet sieving, sedimentation analysis or laser diffraction analysis. Agglomerates are desagglomerated and dispersed in the solution.

The RETSCH ultrasonic baths are also used in chromatography to disperse packing material in the slurry and thus obtain reproducible separation materials.



### **Degassing**

The RETSCH ultrasonic baths are also suitable for degassing solutions or emulsifying oil and aqueous phases.

Order data on page 13


Performance data	UR 1	UR 2	UR 3		
Applications	cleaning, dispersion, degassing				
Feed material	sieves, glass and	d metal compone	nts, suspensions		
Oscillating tank					
ØxH/WxHxD	245 x 130 mm	520 x 200 mm	500 x 300 x 300 mm		
Volumes	5.7 liters	42.0 liters	45.0 liters		
Suitable for the cleaning of	1 sieve	1 sieve	max. 5 sieves		
	200 x 50 mm/8" x 2"	400 x 65 mm	200 x 50 mm/8" x 2"		
Time setting	1 -	15 min or continu	ious		
HF continuous maximum output	2 x 240 W	2 x 600 W	2 x 1000 W		
Technical data	UR 1	UR 2	UR 3		
ØxH/WxHxD	260 x 260 mm	570 x 460 mm	540 x 500 x 340 mm		
Net weight	approx. 5 kg	approx. 21 kg	approx. 27.5 kg		
Noise values (DIN EN 610:	12)				
Noise values	61.5 dB(AU)	76.5 dB(AU)	70.0 dB(AU)		

# Technology of UR 1/2/3

A high-frequency generator produces about 35,000 oscillations per second, which are transferred into the cleaning solution and cause it to resonate. The energy density of the sound field is so high that a cavitation effect sets in. Innumerable extremely small vacuum bubbles develop, which collapse in microseconds due to pressure and suction, in other words they implode. The pulses triggered by this remove dirt particles even at the deepest, least accessible places or they re-

sult in homogenization, dispersion and degassing.

The compact housings and the oscillating tanks are made of stainless steel. Through the ball valve drain located on the casing the cleaning solution can be conveniently and safely drained off. In combination with the splash-proof housing, a high degree of operational safety is provided. Below the oscillating tank is fitted a powerful high frequency generator. The broad-beam oscillation system with PZT oscillators vibrates

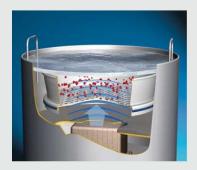

the solvent bath at optimum frequency so as to produce an outstanding cleaning effect.

# Order data DR 100, TG 200, UR 1/2/3

						tem-No.
Vibratory Feed	er DR 100				I	telli-140.
DR 100/40 complete	unit, incl. feeding	kit (push-fit feed chute 40 m	m, holder, hopper 2	.8 liters and fixture)		
DR 100/40 complete	unit for 220-240 V	, 50 Hz			7	0.937.0054
DR 100/40 complete	unit for 110-120 V	, 60 Hz			7	0.937.0055
DR 100 drive unit (p	lease order feeding	kit separately)				
DR 100 drive unit for	220-240 V, 50 Hz				7	0.937.0003
DR 100 drive unit for	110-120 V, 60 Hz				7	0.937.0012
Feeding kits						
Feeding kit with	holder 15/40,	push-fit feed chute 15 mm,	hopper 2.8 liters	and fixture	7	2.020.0001
Feeding kit with	holder 15/40,	push-fit feed chute 40 mm,	hopper 2.8 liters	and fixture	7:	2.020.0002
Feeding kit with	holder 75,	push-fit feed chute 75 mm,	hopper 3.5 liters	and fixture	7:	2.020.0003
Push-fit chutes DR 1	00, lenght 210 mm					
Push-fit feed chute	of stainless steel	, width 15 mm	for holder 15/4	10	0	3.729.0035
Push-fit feed chute	of stainless steel	, width 40 mm	for holder 15/4	10	0	3.729.0036
Push-fit feed chute	of stainless steel	, width 75 mm	for holder 75		0	3.729.0037
Push-fit feed chute	of stainless steel	, width 75/40 mm	for holder 75		0	3.729.0040
Push-fit feed chute	of aluminum,	width 40 mm	for holder 15/4	10	0	3.729.0051
Push-fit feed chute	of aluminum,	width 75 mm	for holder 75		0	3.729.0043
Push-fit feed chute	of stainless steel	, V-shape,	for holder 15/4	10	0	3.729.0039
Holders for push-fit of	thutes DR 100					
Holder 15/40	for push-fit feed	chute 15 mm, 40 mm and V	-shape		0:	3.018.0007
Holder 75	for push-fit feed	chute 75 mm			0	3.018.0008
Hoppers for DR 100						
Hopper 2.8 liters,	of stainless steel	, for chutes 15 and	40 mm		0	3.785.0146
Hopper 3.5 liters,	of stainless steel	for chutes 75/40	and 75 mm		0	2.785.0019
Hopper 0.6 liters,	of stainless steel	for chutes 15 and	40 mm		0	3.785.0151
Hopper 0.6 liters,	of aluminum,	for chutes 15 and	40 mm		0	3.785.0152
	of stainless steel	, for V-chute			0	3.785.0159
Hopper 2.8 liters,						
Fixtures for hoppers	DR 100					
		rs			0	2.266.0167
Fixtures for hoppers	6 liters and 2.8 lite	rs				2.266.0167 2.266.0168
Fixtures for hoppers Fixture for hopper 0.	6 liters and 2.8 lite	rs				
Fixtures for hoppers Fixture for hopper 0.	6 liters and 2.8 lite 5 liters	rs			0.	
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3. Fluid Bed Drye	6 liters and 2.8 lite 5 liters r TG 200	rs id "comfort" with filter bag (	Please order drying	container separately)	0.	2.266.0168
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3. Fluid Bed Drye	6 liters and 2.8 lite 5 liters r TG 200 00, incl. clamping		Please order drying	container separately)	0	2.266.0168
Fixtures for hoppers 0. Fixture for hopper 3. Fluid Bed Drye Fluid Bed Drye TG 2	6 liters and 2.8 lite 5 liters r TG 200 00, incl. clamping V, 50/60 Hz		Please order drying	container separately)	0	2.266.0168 tem-No.
Fixtures for hoppers 6. Fixture for hopper 3. Fluid Bed Drye Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g	6 liters and 2.8 liters  F TG 200  100, incl. clamping IV, 50/60 Hz 200 plass,	id "comfort" with filter bag ( 6 liters	Please order drying	container separately)	0. I 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001
Fixtures for hoppers 5. Fixture for hopper 0. Fixture for hopper 3. Fluid Bed Drye Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of s	6 liters and 2.8 liters  r TG 200  100, incl. clamping V, 50/60 Hz 200 plass, ttainless steel,	id "comfort" with filter bag ( 6 liters 6 liters	Please order drying	container separately)	0 1 7 7 7	2.266.0168 item-No. 0.760.0001 2.783.0001 2.783.0002
Fixtures for hoppers 5. Fixture for hopper 3. Fluid Bed Drye Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g	6 liters and 2.8 liters  r TG 200  100, incl. clamping V, 50/60 Hz 200 plass, ttainless steel,	id "comfort" with filter bag ( 6 liters	Please order drying	container separately)	0 1 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Accessories TG 200	6 liters and 2.8 liters  F TG 200  100, incl. clamping V, 50/60 Hz 200 glass, stainless steel, glass,	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder)		container separately)	0 1 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor	6 liters and 2.8 liters  7 TG 200  100, incl. clamping V, 50/60 Hz 200 glass, stainless steel, glass, t" with filter insert	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep		container separately)	0 1 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of s Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamp	6 liters and 2.8 liters  7 TG 200  100, incl. clamping  V, 50/60 Hz  200  plass,  trainless steel,  plass,  trainless steel,  plass,  trainless steel,  plass,	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep		container separately)	0 1 7 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamp Clamping lid "comfor	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 plass, stainless steel, plass,  t" with filter insert bing lid "comfort", 3 t" with filter bag	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep		container separately)	0 1 7 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamp Clamping lid "comfor Filter bag for clampin	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 plass, stainless steel, plass,  t" with filter insert bing lid "comfort", 3 t" with filter bag ng lid "comfort"	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep		container separately)	7/ 7/ 7. 7. 7. 7. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clampi Clamping lid "comfor Filter bag for clampir Cover with filter inser	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 glass, stainless steel, glass, t" with filter insert bing lid "comfort", 3 t" with filter bag ng lid "comfort" rt for 0.3 liters dry	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces		container separately)	7. 7. 7. 7. 7. 7. 7. 7. 9.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clampi Clamping lid "comfor Filter bag for clampir Cover with filter inser Filter insert for 0.3 li	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 glass, stainless steel, glass, t" with filter insert bing lid "comfort", t" with filter bag ng lid "comfort" rt for 0.3 liters dry ters drying contain	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece		container separately)	7. 7. 7. 7. 7. 7. 7. 9. 0. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clampi Clamping lid "comfor Filter bag for clampir Cover with filter inser Filter insert for 0.3 li Spare drying contain	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 plass, trainless steel, plass, t	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece		container separately)	7. 7. 7. 7. 7. 7. 7. 9. 0. 0.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clampi Clamping lid "comfor Filter bag for clampir Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 10, 50/60 Hz 100 101 101 101 101 101 101 101 101 10	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece		container separately)	7. 7. 7. 7. 7. 0. 0. 0. 0. 0. 0. 0. 0. 0. 0. 0. 0. 0.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping Clamping lid "comfor Filter insert for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 1018ss,	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece		container separately)	7. 7. 7. 7. 7. 7. 7. 0. 0. 0. 0. 0.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 1018ss,  1018 trainless steel, 1018 plass,  1018 trainless steel, 1018 trainless ste	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece		container separately)	7. 7. 7. 7. 0. 0. 0. 0. 0. 0. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping Clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower Quick-clamping elem	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 101ass,	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece		container separately)	7. 7. 7. 7. 0. 0. 0. 0. 0. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower	6 liters and 2.8 liters  T G 200  100, incl. clamping V, 50/60 Hz 200 101ass,	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece		container separately)	7. 7. 7. 7. 0. 0. 0. 0. 0. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower Quick-clamping elem Rods, smooth, 1 pair	6 liters and 2.8 liters  7 TG 200  100, incl. clamping  V, 50/60 Hz  200  glass, stainless steel, glass, tt" with filter insert bing lid "comfort",  "t" with filter bag ng lid "comfort" rt for 0.3 liters dry ters drying containe er of glass, 0.3 liter container 0.3 liters eves 203 mm Ø  10 pieces ents for TG 200, 1	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece		container separately)	7. 7. 7. 7. 7. 7. 7. 0. 0. 0. 0. 0. 0. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of go Drying container of go Accessories TG 200 Clamping lid "comfor Filter insert for clamping Clamping lid "comfor Filter insert for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower Quick-clamping elem Rods, smooth, 1 pair	6 liters and 2.8 liters  T TG 200  100, incl. clamping  V, 50/60 Hz  200  plass,  trainless steel,  plass,  trainless stee	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece pair		container separately)	7. 7. 7. 7. 7. 7. 7. 0. 0. 0. 0. 0. 0. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clampi Clamping lid "comfor Filter bag for clampir Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower Quick-clamping elem Rods, smooth, 1 pair	6 liters and 2.8 liters  T TG 200  100, incl. clamping V, 50/60 Hz  200  plass, plass, trainless steel, plass, trainless steel	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece pair  2 / UR 3 d basket separately)	lacement filters	container separately)	7. 7. 7. 7. 0. 0. 0. 0. 0. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower Quick-clamping elem Rods, smooth, 1 pair  Ultrasonic Batt Ultrasonic baths (ple UR 1 for 230 V, 50/6	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 10, 50/60 Hz 100  101  101  102  103  104  105  105  107  107  107  107  107  107	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm	lacement filters  m, 5.7 liters	container separately)	7. 7. 7. 7. 0. 0. 0. 0. 0. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7.	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying si Dust filter for blower, Quick-clamping elem Rods, smooth, 1 pair  Ultrasonic Batt Ultrasonic baths (ple UR 1 for 230 V, 50/6 UR 1 for 110 V, 60 H	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 17, 50/60 Hz 1800  191ass, 18tainless steel, 18tainless	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 24.5 cm Ø x 13.0 cm	m, 5.7 liters m, 5.7 liters	container separately)	0 1 7 7 7 7 7 7 9 0 0 0 0 0 0 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0001 2.783.0001 2.143.0001 2.143.0001 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001 tem-No. 0.791.0001
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying si Dust filter for blower, Quick-clamping elem Rods, smooth, 1 pair  Ultrasonic Batt Ultrasonic baths (ple UR 1 for 230 V, 50/6 UR 1 for 110 V, 60 H UR 2 for 230 V, 50/6	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 17, 50/60 Hz 1800  1818 glass, 1818 steel, 1818 glass, 1818 it with filter insert 1819 it comfort", 1819 it comfort", 1819 it comfort" 1819 it comfort" 1819 it comfort" 1819 it comfort" 1819 it comfort 1819 it comfo	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece , 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 24.5 cm Ø x 13.0 cm on tank: 52.0 cm Ø x 20.0 cm	m, 5.7 liters m, 5.7 liters m, 5.7 liters m, 42.0 liters	container separately)	0 1 7 7 7 7 7 7 9 0 0 0 0 0 0 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0001 2.783.0001 2.143.0001 2.143.0001 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001 tem-No. 0.791.0001 0.791.0002 0.791.0003
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower, Quick-clamping elem Rods, smooth, 1 pair  Ultrasonic Batt Ultrasonic baths (ple UR 1 for 230 V, 50/6 UR 1 for 110 V, 60 H UR 2 for 230 V, 50 H UR 2 for 110 V, 60 H	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 17, 50/60 Hz 1800  191ass, 18tainless steel, 18tainless	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece rs, 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 24.5 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm	m, 5.7 liters m, 5.7 liters m, 5.7 liters m, 42.0 liters m, 42.0 liters	container separately)	0 1 7 7 7 7 7 7 9 0 0 0 0 0 0 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0001 2.783.0001 2.143.0001 2.143.0001 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001 0.791.0001 0.791.0001 0.791.0003 0.791.0004
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of gone of the second of the seco	6 liters and 2.8 liters  7 TG 200  100, incl. clamping V, 50/60 Hz  200 glass, stainless steel, glass, tt" with filter insert bing lid "comfort", tt" with filter bag ng lid "comfort" ret for 0.3 liters dry ters drying container of glass, 0.3 liter econtainer 0.3 liters eves 203 mm Ø to 10 pieces ents for TG 200, 1  11 SUR 1 / UR tase order cover and to 12 y oscillation to 20, oscillation to 30, oscillation to 40, oscillation to	6 liters 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece rs, 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 50.0 x 30.0 x 30.0 cm on tank: 50.0 x 30.0 x 30.0 x 30.0 cm	m, 5.7 liters m, 5.7 liters m, 5.7 liters m, 42.0 liters m, 42.0 liters cm, 45.0 liters	container separately)	0 1 7 7 7 7 7 7 9 0 0 0 0 0 0 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001 tem-No. 0.791.0001 0.791.0002 0.791.0003 0.791.0004
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of gone of the second of the seco	6 liters and 2.8 liters  7 TG 200  100, incl. clamping V, 50/60 Hz  200 glass, stainless steel, glass, tt" with filter insert bing lid "comfort", tt" with filter bag ng lid "comfort" ret for 0.3 liters dry ters drying container of glass, 0.3 liter econtainer 0.3 liters eves 203 mm Ø to 10 pieces ents for TG 200, 1  11 SUR 1 / UR tase order cover and to 12 y oscillation to 20, oscillation to 30, oscillation to 40, oscillation to	id "comfort" with filter bag ( 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece rs, 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 24.5 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm	m, 5.7 liters m, 5.7 liters m, 5.7 liters m, 42.0 liters m, 42.0 liters cm, 45.0 liters	container separately)	0 1 7 7 7 7 7 7 7 9 0 0 0 0 0 0 7 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.011.0005 2.143.0003 2.737.0003 2.742.0001 0.791.0001 0.791.0002 0.791.0003 0.791.0004 0.791.0005
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 3.  Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container TG Drying container of gone of the second of the seco	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 17, 50/60 Hz 1800  191ass, 18tainless steel, 18tainless	6 liters 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece rs, 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 50.0 x 30.0 x 30.0 cm on tank: 50.0 x 30.0 x 30.0 x 30.0 cm	m, 5.7 liters m, 5.7 liters m, 5.7 liters m, 42.0 liters m, 42.0 liters cm, 45.0 liters	container separately)	0 1 7 7 7 7 7 7 7 9 0 0 0 0 0 0 7 7 7 7 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.011.0005 2.143.0003 2.737.0003 2.742.0001 0.791.0001 0.791.0002 0.791.0003 0.791.0004 0.791.0005
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 1. Fixture for hopper 2. Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying Adapter for drying si Dust filter for blower, Quick-clamping elem Rods, smooth, 1 pair Ultrasonic Batt Ultrasonic baths (ple UR 1 for 230 V, 50/6 UR 1 for 110 V, 60 H UR 2 for 230 V, 50 H UR 2 for 110 V, 60 H UR 3 for 220-240 V, UR 3 for 110-130 V,	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 17, 50/60 Hz 1800  191ass, 18tainless steel, 18tainless	6 liters 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece rs, 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 50.0 x 30.0 x 30.0 cm on tank: 50.0 x 30.0 x 30.0 x 30.0 cm	m, 5.7 liters m, 5.7 liters m, 5.7 liters m, 42.0 liters m, 42.0 liters cm, 45.0 liters		0 1 7 7 7 7 7 7 9 0 0 0 0 0 0 7 7 7 7 7 7 7	2.266.0168  tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.7742.0001 tem-No. 0.791.0001 0.791.0002 0.791.0003 0.791.0004 0.791.0006
Fixtures for hoppers Fixture for hopper 0. Fixture for hopper 1. Fixture for hopper 2. Fluid Bed Dryer Fluid Bed Dryer TG 2 TG 200 for 200-240 Drying container TG Drying container of g Drying container of g Accessories TG 200 Clamping lid "comfor Filter insert for clamping lid "comfor Filter bag for clamping Cover with filter inser Filter insert for 0.3 li Spare drying contain Filter bag for drying si Dust filter for blower, Quick-clamping elem Rods, smooth, 1 pair Ultrasonic Batt Ultrasonic baths (ple UR 1 for 230 V, 50/6 UR 1 for 110 V, 60 H UR 2 for 230 V, 50 H UR 2 for 110 V, 60 H UR 3 for 220-240 V, UR 3 for 110-130 V, Accessories for ultras	6 liters and 2.8 liters  7 TG 200  100, incl. clamping 17, 50/60 Hz 200  18 glass, 18 stainless steel, 18 glass, 18 stainless steel, 19 stainless steel, 10 stainless dry 10 st	6 liters 6 liters 6 liters 3 x 0.3 liters (incl. holder) (exchangeable), incl. 10 rep 10 pieces ing containers, 3 pieces er, 1 piece rs, 1 piece rs, 1 piece pair  2 / UR 3 d basket separately) on tank: 24.5 cm Ø x 13.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 52.0 cm Ø x 20.0 cm on tank: 50.0 x 30.0 x 30.0 cm on tank: 50.0 x 30.0 x 30.0 x 30.0 cm	m, 5.7 liters m, 5.7 liters m, 5.7 liters m, 42.0 liters m, 42.0 liters cm, 45.0 liters cm, 45.0 liters	0249 UR 2: 09.107.0	7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7. 7	2.266.0168 tem-No. 0.760.0001 2.783.0001 2.783.0002 2.002.0005 2.643.0001 2.143.0001 2.107.0002 2.186.0015 2.107.0001 3.186.0024 2.045.0020 2.186.0004 2.001.0005 2.143.0003 2.737.0003 2.742.0001

# **Pellet Press PP 40**



# Benefits at a glance

- Individual pressure force regulation, 1 40 t
- 32 parameter combinations can be stored
- Pressing tools for various diameters and aluminum cups
- Suitable for very hard materials
- Pressure plate made from tungsten carbide
- Easy and safe operation
- Conforming to CE requirements
- 2-year warranty

# Automatic pelletizing for efficient sample preparation


Solid, high-quality pellets are an important precondition for reliable and meaningful XRF analysis. With the PP 40, RETSCH offers a pellet press which presses a wide range of materials such as slag, ores, minerals and cement to strong pellets with a smooth surface. The PP 40 features an **individual pressure force regulation in the range of 1 to 40 t**. Besides controlling the pressure force, it also determines build-up, holding and release of force during pressing. This reduces the inner tensions of the sample and ensures that **even difficult materials are pressed perfectly**.

Performance data	PP 40
Applications	production of pellets for spectral analyses
Feed material	minerals, slag, ores, cement, raw material etc.
Steel rings (external Ø/internal Ø)	51.5 mm/35 mm
	40 mm/35 mm (max. pressure force 20 t)
	40 mm/32 mm
Aluminum cup (external Ø)	40 mm
Parameter combinations	32
Pressure force	1 - 40 t (10 - 400 kN)
Pressure force build-up/holding/release	time respectively 5 - 600 Sek.
Technical data	
WxHxD	836 x 1220 x 780 mm
Net weight	approx. 345 kg
Noise values (Noise measuremen	nt according to DIN 45635-31-01-KL3)
Emission value with regard to workplac	e L <sub>pAeq</sub> 50 dB(A)
	·

## **Technology of PP 40**

The steel ring or aluminum cup is inserted in the pressing tool of the PP 40 and filled with the sample material via a hopper. The steel ring is then pushed beneath the pressure plate and the pressing is started. During pressure build-up the density of the powder increases. The pressure build-up in the PP 40 can be adjusted in such a way that the air inside the hollows of the original powder is pressed out which increases the stability of the pellet. The maximum pressure force must be held over a certain

period of time to allow full development of the interparticulate adhesive forces thus guaranteeing maximum stability. The PP 40 provides pre-selection of the pressure holding time over a period of 600 seconds. During the pressing process the axial movement of the particles at the steel ring produces friction which in turn leads to the formation of a multi-axial stress condition. Therefore, it is important to decrease the pressure evenly and steadily as an abrupt release could lead to the destruction of the pellet.

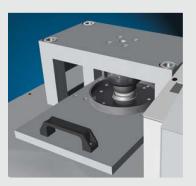

The pressure release time of the PP 40 can be set in such a way that a uniform release of tension is ensured.

### Easy and safe operation

Operation of the PP 40 is very convenient. The pressing tool is mounted on a sliding tray which is pulled out for filling in the sample. The steel ring or aluminum cup is placed into the pressing tool and the material is filled in via a hopper. An advantage of steel rings is lending additional stability to the sample, which is favorable when transporting the sample to the XRF analyzer or in automated systems. Then the pressing tool is pushed beneath the pressure plate made of tungsten carbide, the cover is closed and the pelletizing process is started after selection of the program. The parameters are set easily and safely with one single button, the settings are shown in a graphic display. Up to 32 parameter combinations can be stored with



the PP 40 which guarantees **repro- ducible pelletizing**. When the pressing process is finished, the cover unlocks automatically. The operator can now lift it and pull the pressing tool out to remove the pellet.

The soundproof and completely enclosed Pellet Press PP 40 meets the highest safety standards.

The Pellet Press PP 40 is delivered with mounted pressing tool, of which there are 4 versions:

- for steel rings 51,5 x 8,5 mm, internal diamter 35 mm
- for steel rings 40 x 14 mm, internal diameter 32 mm
- for steel rings 40 x 14 mm, internal diameter 35 mm (max. pressure force 20 t)
- for aluminum cups 40 mm

# **Pellet Press PP 25**



# PP 25 - the "small" solution for XRF analysis

The manual hydraulic Pellet Press PP 25 is a compact benchtop unit with particularly simple and safe operation. With a pressure force of 25 t it is ideally suited for the preparation of solid samples for XRF analysis. The pellets produced are of extremely good quality and are characterized by their high degree of stability. The piston pressure can be read off from the clearly visible manometer scale.

The dies for the Pellet Press PP 25 are available in diameters of 32 mm and 40 mm and can be evacuated completely. This is favorable when pressing porous materials such as e.g. secondary fuels.



## Benefits at a glance

- Produces high quality, stable pellets
- Easy and safe operation
- Pressing tools in 2 sizes; can be evacuated
- Compact benchtop unit
- Conforming to CE requirements
- 2-year warranty

Performance data	PP 25
Applications	production of pellets for spectral analyses
Feed material	minerals, slag, ores, cement, raw material etc.
Dies	32 mm ; 40 mm
Max. pressure force	25 t (250 kN)
Technical Data	
WxHxD	290 x 560 x 310 mm
Net weight	50.5 kg

# Order data PP 40, PP 25

Pellet Press PP 40	Item-No.
Pellet Press PP 40*, complete with die, incl. 5 steel rings or 20 aluminum cups	
Pellet Press PP 40 for 220-230 V, 50/60 Hz, for steel rings Ø 40 /32 mm	20.750.0002
Pellet Press PP 40 for 220-230 V, 50/60 Hz, for steel rings Ø 40 /35 mm	20.750.0003
Pellet Press PP 40 for 220-230 V, 50/60 Hz, for steel rings Ø 51.5/35 mm	20.750.0004
Pellet Press PP 40 for 220-230 V, 50/60 Hz, for aluminum cups Ø 40 mm	20.750.0005
Accessories PP 40	
Steel ring 40 mm outer Ø, 32 mm inner Ø, 1 piece	22.458.0003
Steel ring 40 mm outer Ø, 35 mm inner Ø, 1 piece	22.458.0004
Steel ring 51.5 mm outer Ø, 35 mm inner Ø, 1 piece	22.458.0005
Aluminum cups, straight walls, for tablets with 40 mm diameter, 1000 pieces	22.458.0006
Licowax® C micropowder binder (formerly known as Hoechst Wax), 250 g	22.440.0001
Multimix binder, 500 x 0.25 g tablets	22.440.0002
*Other voltages upon request	

Pellet Press PP 25	Item-No.
Pellet Press PP 25, hydraulic, manual	
(Please order die separately)	20.750.0001
Evacuable dies for Pellet Press PP 25	
Die for 32 mm dia. pellets	22.458.0001
Die for 40 mm dia. pellets	22.458.0002
Accessories PP 25	
Aluminum beaker, sloping walls, for 32 mm dia. pellets, 1000 pcs	22.005.0001
Aluminum beaker, sloping walls, for 40 mm dia. pellets, 600 pcs	22.005.0002
Licowax® C micropowder binder (formerly known as Hoechst Wax), 250 g	22.440.0001
Multimix binder, 500 x 0.25 g tablets	22.440.0002



### Retsch GmbH

Rheinische Straße 36 42781 Haan, Germany

Telephone +49 21 29/55 61 - 0 Telefax +49 21 29/87 02

E-mail info@retsch.com Internet www.retsch.com

a VERDER company

Distributed by:



ADVANCED APPLIED TECHNOLOGIES Contact Us: Irl Ph: 01 4523432

UK Ph: 08452 30 40 30 Web: www.carlstuart.com Email: info@carlstuart.com RETSCH – Your specialist for sample preparation offers you a comprehensive range of equipment. Please request information on our crushers, mills, sieve shakers, sample dividers, feeders as well as cleaning and drying machines.